

ColoQuick Pasteur + MAXX4

HOW TO UPDATE THE SOFTWARE

1. Move the file called "Pasteur.hex" to an empty USB flash drive

2. Insert the USB flash drive

- Pasteur: Left side of the front lid
- MAXX4: Just below the main switch

3. Turn off the machine by using the red main switch

- Wait for the machine to turn off

4. Hold down the "PRG" button while turning the machine back on.



5. Wait for this screen to appear:

- The software is now being updated

```
New application  
Transport init.  
App img. file found_
```

6. The update is completed after 15-30 seconds

```
SW Upgraded
```

7. This screen will appear after a few seconds, showing the updated software version.

- In this example, version 2.47a.

```
Pasteur Maxx 4  
ver. 2.47a / 251  
BSL: 5
```

The ColoQuick unit is now ready for use, and the USB flash drive can be removed.

